

SpinningDisc Switch On/Off Procedures

1. Switch on Microscope (wall plug labelled Power Strip 1 of 3)
2. Switch on Spinning Disc Boxes (wall plug labelled Power Strip 2 of 3)
3. Switch on Lasers (only when you use lasers, wall plug located behind laser box labelled Lasers & Confocal)
4. Turn the laser safety key to ON (labelled Safety Shutter on laser boxes)
5. Switch on camera (at the back panel of the camera)
6. Log in
7. Run SlideBook6

If you need to do TIRF experiment, please ask Yan in advance.

Need help? Contact Yan Gu: 07981875417, office: 2c37, ext:6997

1. Close SlideBook
2. Log off
3. Switch off camera (on back panel of the camera)
4. Turn the laser safety key to OFF (labelled Safety Shutter on laser boxes)
5. Switch off lasers (only when switched on, wall plug labelled Lasers & Confocal)
6. Switch off Spinning Disc Boxes (wall plug labelled Power Strip 2 of 3)
7. Switch off microscope (wall plug labelled Power Strip 1 of 3)
8. Clean the objective (with lens tissue)