

Apotome Switch On/Off Procedures

1. Switch on Heater & CO₂ (only for live imaging, wall plugs labelled 'Heater' & 'CO2')
2. Switch on Apotome (wall plug labelled 'Apotome')
3. Switch on Microscope (back-left side panel on microscope frame, labelled 'Microscope')
4. Switch on CameraT (wall plug labelled 'CameraT')
5. Switch on CameraR (only for live imaging, wall plug labelled 'CameraR')
6. Log in & run Zen (icon with blue background)

1. Exit Zen & log off
2. Switch off CameraR (only if it is on, wall plug labelled 'CameraR')
3. Switch off CameraT (wall plug labelled 'CameraT')
4. Switch off Apotome (wall plug labelled 'Apotome')
5. Switch off Heater & CO₂ (only if it is on, wall plugs labelled 'Heater' & 'CO2'))
6. Clean the objective (with lens tissue)

If you need to do dual camera live experiment, please ask Yan in advance

Need help? Contact Yan Gu: 07981875417, office: 2c37, ext:6997