

AxioVert Switch On/Off Procedures

1. Check that the computer is ON.
2. Switch on microscope (Wall plug labelled [Stripe 3](#))
3. Switch on camera (camera controller labelled [Camera on shelf](#))
4. Switch on fluorescence lamp (HXP120V lamp box on shelf labelled [Lamp](#))
5. Log into your university account
6. Run Micro-manager

1. Close Micro-manager
2. Log off
3. Switch off fluorescence lamp (HXP120V lamp box on shelf labelled [Lamp](#))
4. Switch off camera (camera controller labelled [Camera on shelf](#))
5. Switch off microscope (Wall plug labelled [Stripe 3](#))
6. Remove your slides and clean the objective with lens tissue

Need help? Contact Yan Gu: 07981875417, office: 2c37, ext:6997