

# Dragonfly Switch On/Off Procedures

1. Switch on heater and 5% CO<sub>2</sub> (only when imaging live samples)
2. Check Workstation is on
3. Switch on Stripe1 (labelled on wall plug)
4. Switch on Stripe2 (labelled on wall plug)
5. Switch on Stripe3 (labelled on wall plug)
6. Login university account
7. Run Fusion software

1. Close Fusion & log off
2. Switch off Stripe3 (labelled on wall plug)
3. Switch off Stripe2 (labelled on wall plug)
4. Switch off Stripe1 (labelled on wall plug)
5. Turn off heater and 5% CO<sub>2</sub> (only when imaging live samples)
6. Clean the objective (with lens tissue, only applied to water/oil objective)

If you plan live or selective illumination experiment , please ask Yan in advance.

**Need help?** Contact Yan Gu: 07981875417, office: JMS 2c37, ext:6997