

E400 Switch On/Off Procedures

1. Make sure the workstation is on
2. Make sure the camera is on (LED indicator on the back panel of the camera is ON)
3. Switch on fluorescence light source (ON/OFF button on the light control pad)
4. Log into university account
5. Start MicroManager software

1. Close MicroManager software
2. Log off university account
3. Switch off fluorescence light source (wall plug labelled light)
4. Clean the objective if you used Oil objective (use lens tissue only)

Need help? Contact Yan Gu: 07981875417, office: JMS 2c37, ext:6997