

IX70 Switch On/Off Procedures

1. Switch on Stage (wall plug labelled Stage)
2. Switch on fluorescence light source (on the shelf)
3. Switch on CCD camera (on the front panel of camera drive on the shelf)
4. Log in
5. Run Micro-Manager program (configure file (*.cfg) is in C:\temp)

1. Close Micro-Manager Program
2. Log off
3. Switch off CCD camera (on the front panel of camera drive on the shelf)
4. Switch off fluorescence light source (keep it on if anybody booked behind you within 1 hour)
5. Switch off Stage (wall plug labelled stage)
6. Clean the objective (with lens tissue)