

# LSM880 Switch On/Off Procedures

1. Switch on Main Switch (labelled '1' on switch pad)
2. Switch on Systems/PC (labelled '2' on switch pad)
3. Switch on Components (labelled '3' on switch pad)
4. Check Laser key is ON (right side on switch pad)
5. Check laser key is at '-' position (front panel of LASOS laser driver)
6. Log in
7. Run Zen (icon with black background)

1. Shut down lasers from software
2. Close Zen & log off
3. Switch off Components (labelled '3' on switch pad)
4. Switch off Systems/PC (labelled '2' on switch pad)
5. Wait for the cooling fans to stop (noticeable reduced noise level from LASOS laser driver)
6. Switch off Main Switch (labelled '1' on switch pad)
7. Clean the objective (with lens tissue)

If you need to do live experiment, please ask Yan in advance.

**Need help?** Contact Yan Gu: 07981875417, office: 2c37, ext:6997