

# Operetta Switch On/Off Procedures

1. Make sure the workstation is on
2. Switch on the power of Operetta  
(Green switch on the back of left side of microscope)
3. Log into university account
4. Wait until Operetta's initialization finishes
5. Start **Harmony** software

## If it is a live experiment:

- Switch the heater on from **Harmony** 3 hrs in advance
- Turn on CO<sub>2</sub> before experiment

1. Take out the sample
2. Press **Load** button in **Harmony** to close the sample loading chamber
3. Close **Harmony** software
4. Log off university account
5. Switch off **the power of Operetta**  
(Green switch on the back of left side of microscope)

## If it is a live experiment:

- Turn off CO<sub>2</sub>