

SP8 Switch On/Off Procedures

1. Make sure the workstation is on
2. Switch on the **Laser Power** (On the switch panel left hand underneath table)
3. Switch on the **Scanner Power** (On the switch panel left hand underneath table)
4. Switch on **PC Microscope** (On the switch panel left hand underneath table)
5. Turn the **Laser Emission** key to **On-1** (On the switch panel left hand underneath table)
6. Switch on fluorescence lamp (ET6000 on shelf)
7. Log into university account
8. Start **LAS X** software

1. Close **LAS X** software
2. Log off university account
3. Switch off **PC Microscope** (On the switch panel left hand underneath table)
4. Switch off **Scanner Power** (On the switch panel left hand underneath table)
5. Turn off **Laser Emission** key to **Off-0** (On the switch panel left hand underneath table)
6. Wait for the cooling fan to stop (noticeable reduced noise from laser fan)
7. Switch off Laser Power (On the switch panel left hand underneath table)
8. Clean the objective (with lens tissue, only applied to oil objective)