

# Abberior STED Switch On/Off Procedures

To achieve good resolution, please switch on the system 2 – 3 hours in advance

1. Check the workstation is ON
2. Switch on Microscope (wall plug labelled Microscope)
3. Switch on fluorescence light source (wall plug labelled Light)
4. Switch on Microscope Control (on the back panel of Control Hub)
5. Log in
6. When initialization of Control Hub finishes, run *Inspector* software

1. Close *Inspector* software
2. Log off
3. Switch off fluorescence light (wall plug labelled Light)
4. Switch off Microscope (wall plug labelled Microscope)
5. Remove sample and clean the objective (with lens tissue)

**Need help?** Contact Yan Gu: 07981875417, office: JMS 2c37, ext:6997  
or Office: CRPC 3.15, ext 3908